

Rodent Warmer

by Stoelting

Cat. No. 21100-800

General

Use Rodent Warmer before, during and after surgical procedures to improve surgical outcome and overall longevity.


Monitor and maintain animal temperature with three programmable settings: animal specific, timer or use with thermal probe.

The Rodent Warmer is available with mouse, rat or home-cage heating pad; the rectal thermal probe is sold separately.

The Rodent Warmer is a perfect complement to our line of Ventilators and Anesthesia Systems, as well as our BP Recorder, and Stoelting's Stereotaxic Instruments.


Model X2 is also available, which allows for easy programming and controls 2 independent heating pads simultaneously.

Available with a combination of mouse, rat or home-cage heating pads.

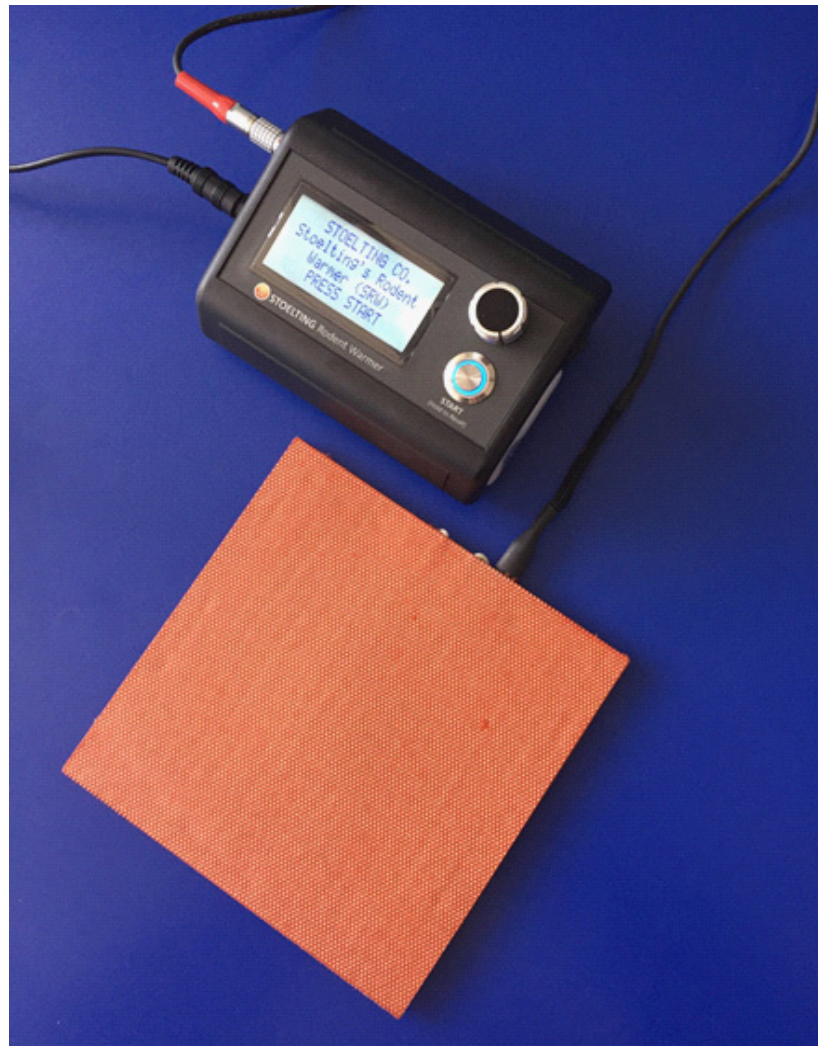

Minimize heat loss and improve surgical outcome with the new rodent warmer!

**Compact
for Mice and Rats**

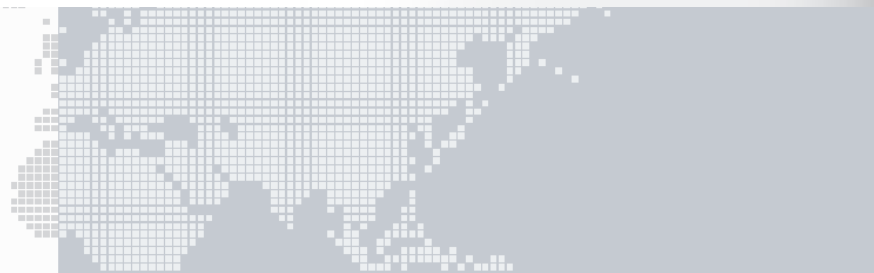

Main Features

- Ideal for use with mice and rats
- Easy to use dial control
- Heating Pad Included (mouse, rat, or home cage)
- Preprogrammed Animal Temperatures
- Lightweight, small footprint
- Use with or without rectal probe (*rectal probe sold separately*)

General

The new Rodent Warmer can be used as a general warming system or with a rectal probe (sold separately) for more accurate, core temperature monitoring during pre- and post-op surgical procedures.

Before: Heating pad can be placed underneath an induction chamber to reduce heat loss during anesthetic administration.

During: Place heating pad on a rodent surgery table or stereotaxic instrument to maintain and monitor temperature during surgical procedures, ventilation and anesthesia.

After: Cage heating pads can be placed in the animal's home cage for faster recovery following surgical procedures.

Instrument Description

The Rodent Warmer provides three Operating Settings:

1. Select from pre-programmed animal temperatures
2. Countdown Timer
3. Use with Thermal Rectal Probe

All selectable via an easy-to-use dial control.



Heating Pads

Heating pads are available in three different sizes:



Cage Heating Pad
16x38cm

Rat Heating Pad
15.25x15.25cm

Mouse Heating Pad
7x7cm

Optional

Rodent Warmer X2 is also available, to control 2 independent heating pads simultaneously. Available with a combination of mouse, rat or home-cage heating pads.



Physical

- Temperature Control Range : 25-45°C
- Temperature Resolution : 0.1°C
- Heater Blanket Connection : 4-pin locking DIN
- Heater Power : 24VDC@3A
- Max Power Requirements : 120/240VAC (switchable)
50/60Hz, 75VA
- Probe Input Connector : Phone jack
- Probe Dimensions : 0.62in/1.6mm tip diam.
- Control Box Dimensions : 12.5(l)x9.5(w)x4(h)cm
- Control Box Weight : 200g

Ordering Information

Rodent Warmer X1

- 21100-800M Rodent Warmer with Mouse Heating Pad
- 21100-800R Rodent Warmer with Rat Heating Pad
- 21100-800C Rodent Warmer with Cage Heating Pad

Rodent Warmer X2

- 21100-850MM Rodent Warmer with 2 Mouse Heating Pads
- 21100-850RR Rodent Warmer with 2 Rat Heating Pads
- 21100-850CC Rodent Warmer with 2 Cage Heating Pad

Accessories:

- 21100-810 Mouse Heating Pad
- 21100-812 Rat Heating Pad
- 21100-814 Cage Heating Pad
- 21100-304 Rectal Thermal Probe (*)

(*) Rectal Thermal probe always sold separately.